

Anti- R-FNR (Root Ferredoxin-NADP reductase) antibody, rabbit polyclonal 81-007 100 µg

Shipping and Storage: Shipped at 4°C or -20°C and store at -20°C. Do not freeze.

Immunogen: Purified recombinant maize root-FNR protein (full-size, no-tag attached)

Form: 1 mg/ml in PBS- with 50% glycerol. Filter sterilized. Azide and carrier free.

Purity: IgG, ProteinA-affinity purified from rabbit antiserum

Reactivity: Plant root FNR proteins (R-FNR1 and R-FNR2) including Arabidopsis and Maize. The antibody also reacts with leaf FNRs.

Applications:

- 1. Western blotting (1/1,000~1/30,000 dilution)
- 2, ELISA (assay dependent)

Other applications have not been tested.

Background: Ferredoxin-NADP reductase (FNR) isoproteins of plant roots play a key role in redox homeostasis of NADPH / NADP+ and donation of reducing equivalence to ferredoxin. R-FNR2 is major form of R-FNR and involved in reduction and detoxication of nitrite in root.

Data Link: UniProtKB: Q41736 (Z. mays), Q9M0V6 (A. thaliana)

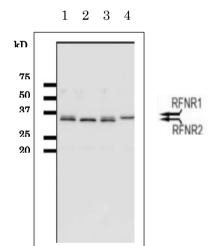


Fig.1 Detection of Arabidopsis R-FNR1 and R-FNR2 isoproteins by western blotting with anti-R-FNR antibody Extracts from wild type strain Arabidopsis Col (1 and 3), mutant rfer1 (2) and rfnr2-2 (4) grown under 0.2 mM nitrate for 7 days were analyzed by western blotting. Anti-R-FNR antibody was used at 1/2,000 dilution. Note that R-FNR2 is dominant form in wild-type roots. Wild type produces both R-FNR1 and R-FNR2. Mutant rfnr1 produces R-FNR2 and mutant rfr2 produces R-FNR1.

Reference: This product has been used in the following publication.

- Onda Y, et al. Differential interaction of maize root ferredoxin:NADP(+) oxidoreductase with photosynthetic and non-photosynthetic ferredoxin isoproteins. Plant Physiol. 2000, Jul;123(3):1037-45. PMID: 10889253 WB; Maize
- Hachiya T et al. Arabidopsis Root-Type Ferredoxin:NADP(H) Oxidoreductase 2 is Involved in Detoxification of Nitrite in Roots. Plant Cell Physiol. 2016 Nov;57(11):2440-2450. PMID: <u>27615794</u> WB; Arabidopsis



BioAcademia,Inc. Tel. 81-6-6877-2335 Fax. 81-6-6877-2336 <u>info@bioacademia.co.jp</u> <u>http://www.bioacademia.co.jp/en/</u>